

岛津SHIMSEN系列 过滤式样品瓶

SHIMSEN Mini-Filter Vial





采用SHIMSEN Mini-Filter vial 过滤样品瓶可代替针式滤器、注射器、样品瓶三个不同的实验组件，可以简化工作流程，减少实验室中的浪费，帮助实验人员提升实验效率。

- 过滤完的样品可以直接上机实验，减少不必要操作，节约时间成本；
- 减少处理过程中与样品接触，样品受污染几率大大减少；
- 不同颜色盖子对应不同类型的膜；




■ 如何使用？

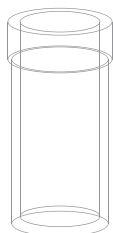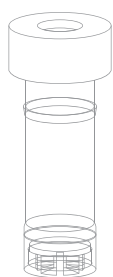


移入样品

用移液器添加样品到
填充线上

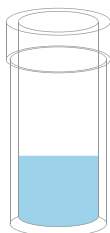

按压

向下轻轻按压

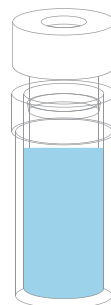

按压至底部

缓缓将带有滤膜的内部小瓶按
压至底部

SHIMSEN Mini-Filter Vial

P01-P03



380-06211-01



380-06211-02



380-06211-03



380-06211-04

产品编号	380-06211-01	380-06211-02	380-06211-03	380-06211-04
产品描述	Nylon 0.22 μm Clear Filter Vial,pp	Nylon 0.45 μm Clear Filter Vial,pp	PTFE 0.22 μm Clear Filter Vial,pp	PTFE 0.45 μm Clear Filter Vial,pp
100 pcs/pk.				



380-06211-05



380-06211-06



380-06211-07



380-06211-08

产品编号	380-06211-05	380-06211-06	380-06211-07	380-06211-08
产品描述	PES 0.22 μm Clear Filter Vial,pp	PES 0.45 μm Clear Filter Vial,pp	PVDF 0.22 μm Clear Filter Vial,pp	PVDF 0.45 μm Clear Filter Vial,pp
100 pcs/pk.				



380-06211-09



380-06211-10



380-06211-11



380-06211-12

产品编号	380-06211-09	380-06211-10	380-06211-11	380-06211-12
产品描述	Nylon 0.22 μm Amber Filter Vial,pp	Nylon 0.45 μm Amber Filter Vial,pp	PTFE 0.22 μm Amber Filter Vial,pp	PTFE 0.45 μm Amber Filter Vial,pp
100 pcs/pk.				

SHIMSEN Mini-Filter Vial

SHIMSEN Mini-Filter Vial

P01-P03



380-06211-13



380-06211-14



380-06211-15



380-06211-16

产品编号	380-06211-13	380-06211-14	380-06211-15	380-06211-16
产品描述	PES 0.22 μm Amber Filter Vial,pp	PES 0.45 μm Amber Filter Vial,pp	PVDF 0.22 μm Amber Filter Vial,pp	PVDF 0.45 μm Amber Filter Vial,pp
100 pcs/pk.				



380-06211-17



380-06211-18



380-06211-19



380-06211-20

产品编号	380-06211-17	380-06211-18	380-06211-19	380-06211-20
产品描述	Ultra-clean PTFE 0.22 μm Clear Filter Vial,pp	Ultra-clean PTFE 0.45 μm Clear Filter Vial,pp	Ultra-clean PTFE 0.22 μm Amber Filter Vial,pp	Ultra-clean PTFE 0.45 μm Amber Filter Vial,pp
100 pcs/pk.				

尺寸：12mm，34mm；

是否灭菌：否

过滤面积：0.28cm²

最大装样量/滞留体积：480ul/30ul

最大压力：1.5 bar (21.6 psi)

最高工作温度：121 °F (50 °C)

标柱颗粒保留效率 $\geq$ 95%



Shimadzu (Shanghai) Global Laboratory Consumables Co.,Ltd.

www.sglc.shimadzu.com.cn

www.shimadzumall.com

Contact: contact@sglc.shimadzu.com.cn

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.